# KAME ver 4.1.10 (Partially in English)

Sep 23, 2015 Kentaro Kitagawa

## Preface

---

“kame” is a C++ multithreaded program which focuses on automated measurement by scripting with Ruby language.

It is especially intended to be used in solid-state NMR experiment, but includes many types of measurement drivers. The source is cross-platform (Linux/Mac OS X/Windows).

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# To install kame on Mac/Linux machines

See japanese document or memos in the source zip.

# To install kame on Windows machines

## The case of using compiled binaries

Qt of version 5.3 or later including MinGW32bit is necessary.

If you need GPIB,

Install a 488.2 driver from National Instruments (support for ANSI C and direct entries are needed)

If you need DAQmx,

Install a DAQmx driver from National Instruments (support for ANSI C is needed)

Expand kame-4\*\*-win32.zip and launch it via kame.bat. Use not too old machine so that it has enough functioning support for OpenGL2.1.

If you use USB devices of Thamway, copy fx2fw.bix, slow\_dat.bin, fullspec\_dat.bin to the directory of kame executable. If something is wrong, try deleting fullspec\_dat.bin.

## The case of compiling source zip

First, try the above instructions.

Next, several depending libraries.

fftw-3.3.4-dll32.zip (Expand it in the upper directory of the kame source, and name the directory "fftw3")

<http://www.fftw.org/download.html>

Install Ruby and msys (<http://www.mingw.org/>), and activate the path for the ruby. And expand ruby-2.2.\* source in the upper directory of the kame source, and change the name to "ruby".

Launch msys.bat, and in the ruby source directory, activate the path of the mingw (suchlike... PATH=\$PATH:/c/QT/Qt-(somewhat)/Tools/mingw4(somewhat)\_32/bin), and type

```
CFLAGS="-g0" ./configure --enable-shared --disable-rubygems
```

```
make
```

Then you would obtain a DLL of Ruby

For zlib, get "Compiled DLL" from the following, and expand it in the upper directory as well and name it "zlib"

<http://www.zlib.net/>

You need to compile the newer version of GSL by yourself

<http://www.gnu.org/software/gsl/>

Download the newest gsl, and expand it in the upper dir, naming "gsl". Do the following as you have done for ruby

```
./configure;make
```

Then, copy ./libs/libgsl-0.dll and cblas/.libs/libgsl-cblas-0.dll to the directory of the executables.

Open kame.pro from Qt creator. Build it in the mingw32 environment.

To execute kame, you need to copy the following dlls to the directory of the kame executables (for example, build-kame-\*\*\*-Debug).

zlib1.dll(in C:\Program Files\*\GnuWin32\bin or so), llibgsl.dll, libgslcblas.dll (in gsl) ,  
ibftw3-3.dll (in FFTW3) , msvc\*-ruby\*.dll(in ruby)

And you can launch kame from Qt Creator.

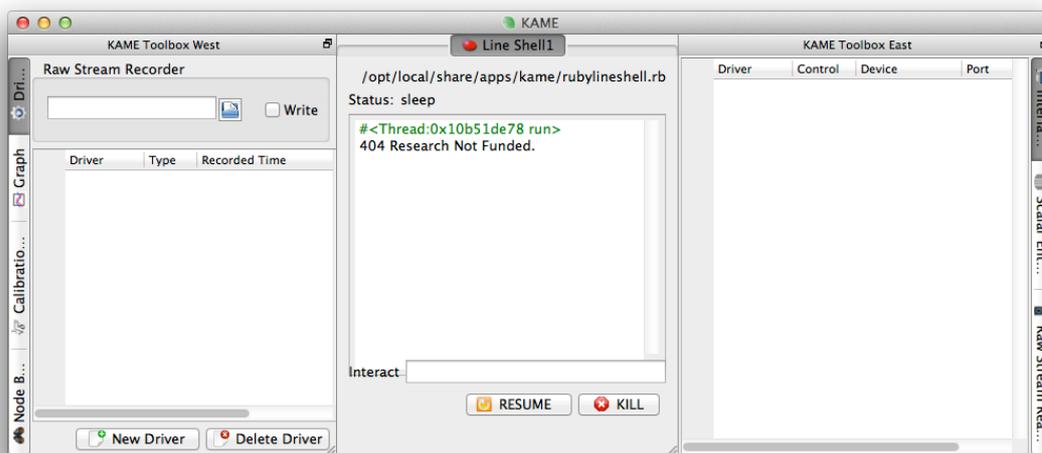
For i18n, kame\_ja.qm must be in the same directory, and “kame/script/rubylineshell.rb” is needed in the “Resources” directory to use a line shell.

## First step

---

Error logs will be written in /tmp/kame.log(Unices) or in the current directory(Windows).

## Basic usage

(First time) Add favorite driver for your instrument by invoking “New” button in the “Driver” tab.

(Next time) Open your .kam file to restore the saved state in the menu “File”-“Open”

Push “Start” button in the “Interface” tab. Before this, you must specify address and/or port for GPIB/SerialPort/TCP.

When you click a label of driver in “Driver” or “Interface” tab, you can open the driver specific window

The tab “Scalar Entry” is intended to prepare the conditions for storing line-by-line text data. The entries would be temperature or voltage etc.. When the difference after the last written value exceeds “Delta”, one line will be appended in the file. If “Delta” is negative, it is appended on getting new value. When you click the cell, a chart as a function of time will appear.

The tab “Graph” is for making X-Y-(Z) graphs from scalar entries.


To save the state in kame, invoke “File”-“Save”. Frequent saving may save your time.

## Usage for text boxes

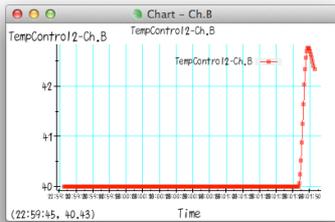When you edit the text of UIs, you have to press “Enter” key at the end, because the instruments should not reflect the intermediate text. Unless so, the original state is recovered.

The values in most of the text boxes can be operated from the Ruby scripts.

## Usage for Graphs/Charts

They are drawn by OpenGL in 3D space as transparent entities.

Left double click opens detailed dialog.

Right double click opens hints for you.

Left click and dragging on the plot can be used to expand scales.

Right click on the plot enables auto scaling.

Wheeling on the plot is zooming in or out.

Right click and dragging on the axis expands the axis.

Right click on the axis enables auto scaling for the axis.

Wheeling on the axis inclines the plot in the 3D space.

Middle click and dragging inclines the plot in the 3D space more freely.

Middle click restores a view point.

If you are dealing with the wave data, which can be stored in the text, you have a filepath bar above or beneath the graph. The directory icon can choose the file name, but you must not forget to click “Dump” when you want to save the data.

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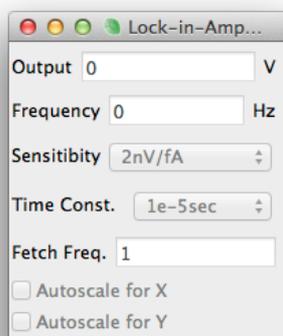
Stanford Resresearch SR830 lock-in amplifier (GPIB)

NF LI5640 lock-in amplifier (GPIB)

Andeen-Hagerling 2500A capacitance bridge (GPIB)

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
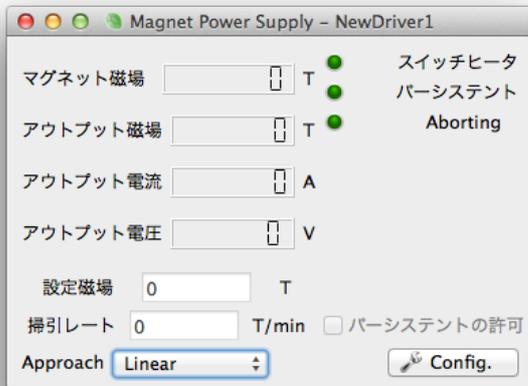
Oxford PS/IPS-120 magnet power supply (GPIB, SerialPort)

Cryogenic SMS10/30/120/150C magnet power supply (USB Serial Port)

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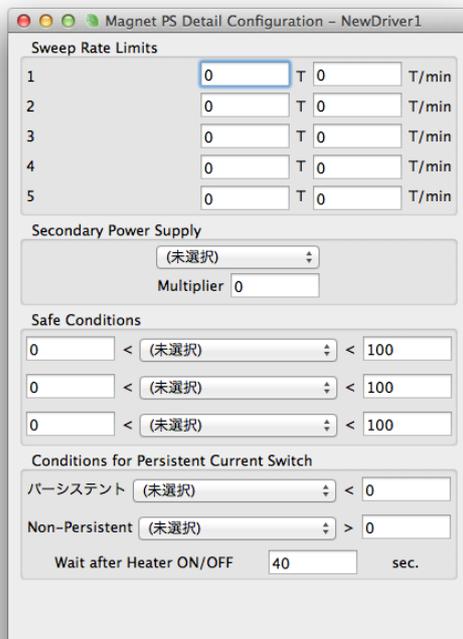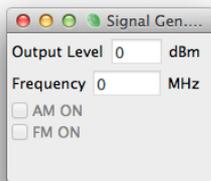
DSTech. DPL-3.2XGF (SerialPort)

HP/Agilent 8643/8644/8648/8664/8665 signal generator (GPIB)

Rhode-Schwartz SML-01/02/03/SMV-03 signal generator (GPIB, SerialPort)

Thamway NMR PROT built-in signal generator (TCP/IP, GPIB, USB)

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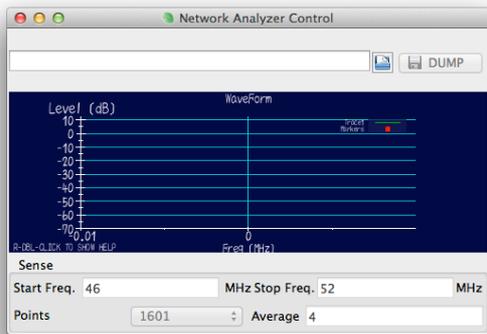
HP/Agilent 8711/8712/8713/8714 network analyzer (GPIB)

Agilent E5061/E5062 network analyzer (GPIB)

DG8SAQ VNWA3E w/ custom dll network analyzer (TCP/IP)

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Cryocon M32/M62 temperature controller (GPIB)

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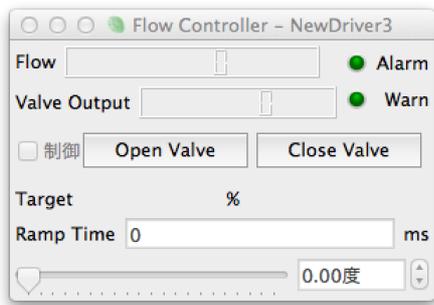
Fujikin FCST1000 Series Mass Flow Controllers (Serial Port 38400bps)

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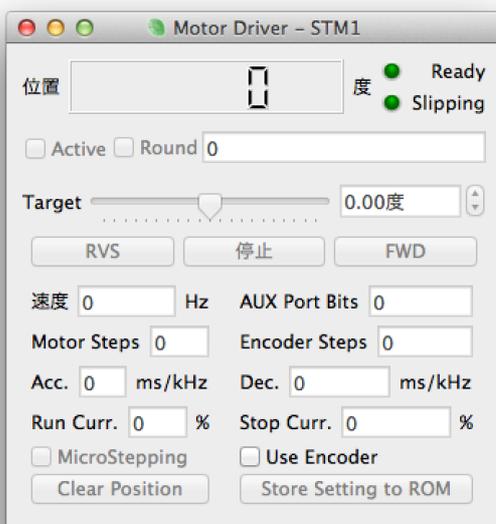
## NMR Pulser

NMR pulser on NI-DAQ M series (NI-DAQmx)

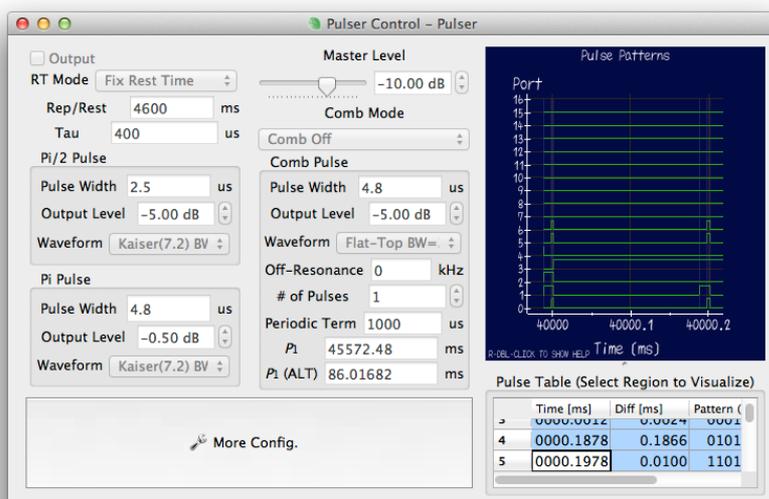
NMR handmade pulser on H8 (SerialPort)

NMR handmade pulser on SH2 (SerialPort)

NMR pulser Thamway N210-1026 PG32U40 (USB)

NMR pulser Thamway N210-1026S/T (GPIB, TCP/IP: experimental)

See appendix for DAQmx

The pulser drivers without featuring QAM cannot change the items of intensities and waveforms.

Digital waves will appear right side.

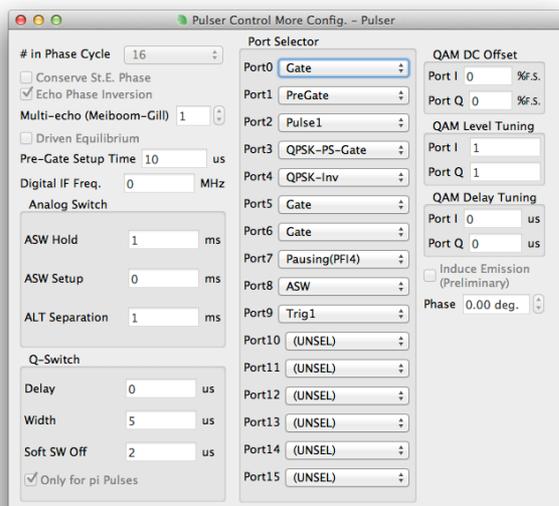
When "RT mode" is "Fix Rep. Time", the total period between the 1st (Pi/2) pulse is fixed to "Rep/Rest" time. Otherwise, ("Fix Rest Time"), the clearance time between the last pulse or  $2\tau$  to Saturation (Comb) Pulse is fixed to the described value.

The times (suchlike "Tau", " $P_1$ ") are defined to the center of pulses, in principle. For "Comb pulse", the center of the last saturation pulse is considered.

When "Comb mode" is "P1 ALT", it uses two  $P_1$  values alternatively. In that case, two FID/echo drivers are necessary to acquire  $T_1$ . Hardware gating using "ASW" signal is necessary.

Seemingly, "Comb ALT" mode alternatively issues the pulse sequence with and without saturation pulses.

The following window appears when "More Config." button is invoked.

“# in Phase Cycle” determines the cycles for  $\pi/2, \pi$  pulses up to the value of 4, and in addition, the phases for comb pulses can be altered when the # > 4.

“Conserve St.E. Phase” allows to measure Stimulated Echo with phase cycling. For this, use two pulses for the comb, and the period of the comb should be the same of  $\tau$ , and width of  $\pi/2$  pulse is zero, and the width of  $\pi$  pulse should be same of that of comb.

“Multi-echo” more than 2 uses Carr-Purcell-Meiboom-Gill (CPMG) cycles. It features phase cycling even for CPMG. For this, you have to do proper setting in the FID/echo driver.

“Port selector” is used for associating the hardware ports and the functions.

“Gate” is gating.

“PreGate” is gating with a setup time for a power amp.

“Trig1” is a trigger rising at  $2\tau$ .

“Trig2” is a trigger rising at  $\pi/2$  pulse and falling at  $\pi$  pulse.

“ASW” is gating for the receiver.

“QSW” becomes H level only after the pulse. This is for Q switch.

“Pulse1” is identical to  $\pi/2$  pulse.

“Pulse2” is identical to  $\pi$  pulse.


See QPSK logics below:

(Phase)	“QPSK-A”	“QPSK-B”	“QPSK-NonInv”	“QPSK-Inv”	“QPSK-PS-Gate”
0°	L	L	H	L	L
90°	H	L	H	L	H
180°	H	H	L	H	L
270°	L	H	L	H	H

Thamway receiver would use a pair of “QPSK-A”, “QPSK-B” or “QPSK-PS-Gate”, “QPSK-Inv”

## NMR FID/Echo measurement

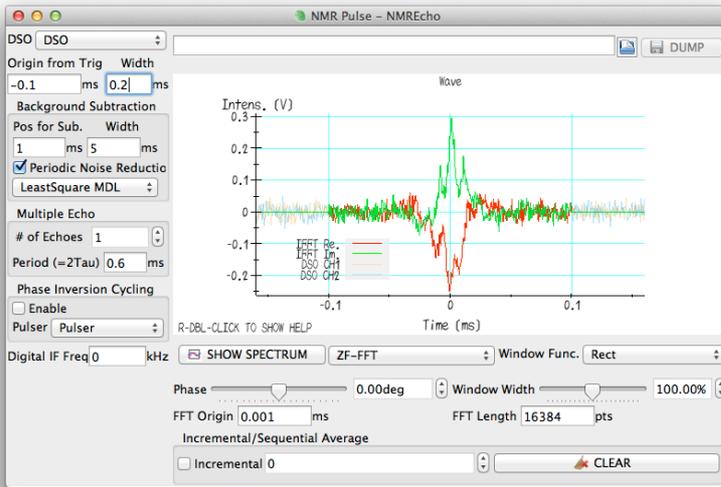This driver is an analyzer, which subtracts a background level from the DSO wave, cuts it using a window, sums up, and makes a frequency-domain spectrum.

For ordinary echo measurement, set the trigger of DSO to  $2\tau$  (= "Trig1"), and use "Single Sequence" mode for DSO.

It cuts the wave of DSO from the "Origin from Trig." by "Width". But for "Multiple Echo", displayed waves are averaged ones for "# of Echoes" times.

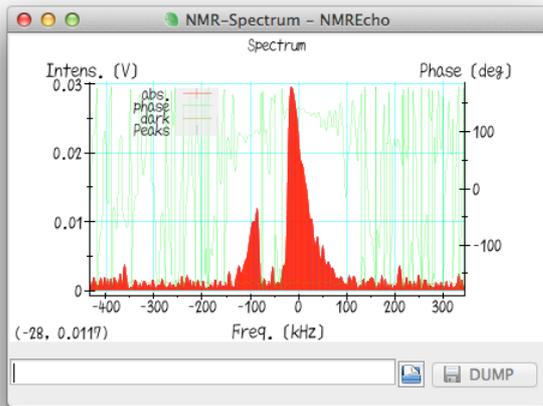
When "Width" of "Background Subtraction" is not zero, it displays a subtracted data using the average level from "Pos for Sub." with "Width". For better S/N, you have to use much larger "Width" than that of "Width" for echoes.

If "Period Noise Reduction" is activated, a least square fitting for periodic noise in the background part is attempted to extrapolate the noises. The number of noise frequencies are estimated by the information criterion. This function requires a lot of CPU powers.

When the "Phase Inversion Cycling" is on, it averages a positive and negative echoes alternatively after an inversion for the latter. This procedure allows to cancel out the time-dependent background.

When the "Incremental" is activated, once the stored wave is cleared and it averages all waves forever unless you invoke "CLEAR" button. The number box displays the average count.

Clicking "SHOW SPECTRUM" puts the following window.

Usually we use "ZF-FFT" (Zero-Filling Fast Fourier Transform) mode. The origin for the FFT is given by "FFT Origin", which is counted from the trigger. Of course, you can use "Window func."


The red bars are the absolute in the complex FFT. The phase (green) should be flat if the "FFT Origin" is proper. "dark" is a noise level which is estimated by the background. "Peaks" are maxima in "abs."

The displayed "IFFT Re. / IFFT Im." in the main window are IFFT result for the spectrum analysis. In ZF-FFT, it is nothing but a wave after windowing. "DSO Re. / DSO Im." is the waves of DSO after the background subtraction.

In the scalar entries, the maximum value and its frequency is emitted.

## NMR Relaxation Rate Measurement

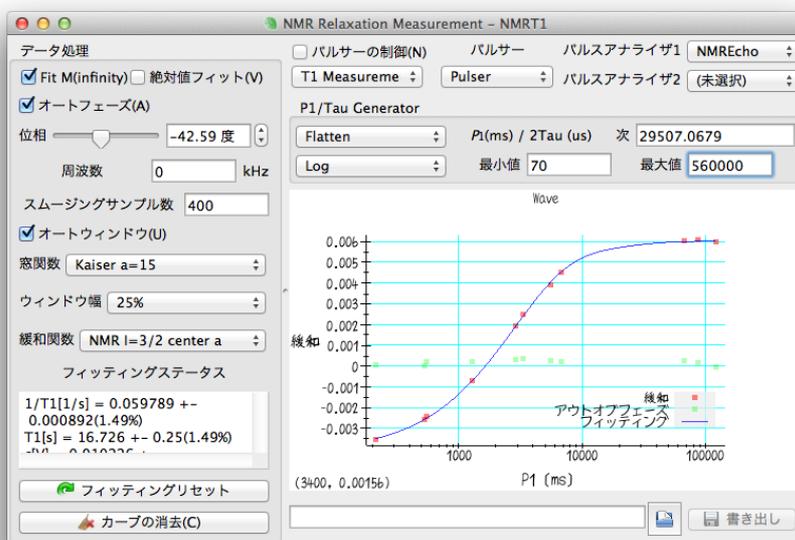The driver is analyzer, which measures relaxation rates using waves in FID/Echo drivers. Of course the pulser is controlled by this. Configurations for window function are not shared with FID/Echo drivers, and can be freely changed during your measurement.

Invoking “Control Pulser” first clears the results and new  $P_1$  value (written in “Next”) will be transferred to the pulser driver when a new data come in. When the generator is “Random”,  $P_1$  will be randomly chosen between “Min” and “Max”. The “Flatten” generator tries to find the most vacant region. “Log” distribution is standard for the generators. You can change “Next” value for  $P_1$  at any time.

“Fitting Status” summaries a result of a least square fitting for a relaxation curve.

“Auto Phase” choses automatically the phase that can maximize the difference in the curve.

“Smooth Samples” specifies a number of regions for the truncation for  $P_1$  before fitting.

“Auto Windows” choses automatically the window function that can maximize S/N. It is not suited for an experiment where a frequency dependence of  $T_1$  is physically important.

“Fit M(infinity)” adds a fitting parameter for a value at  $t \rightarrow \infty$ . This should be turned off for  $T_2$  measurement.

“Frequency” is a center frequency for the relaxation analysis. Change of this entity clears the stored result.


“Reset Fitting” could be used when the fitting is not well functioning due to the initial values.

$1/T_{1,2, \text{st.e.}}$  values and errors will be emitted to the scalar entries.

## NMR Spectrum Measurement (Frequency Sweep)

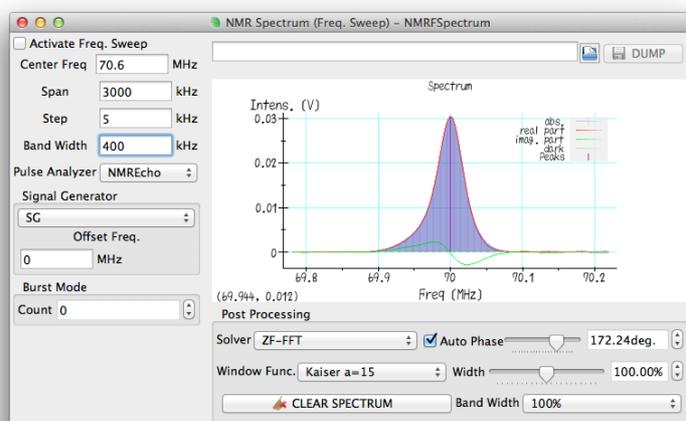This driver is an analyser which measures frequency-swept spectrum using waves in FID/Echo driver. Configurations for window function are not shared with FID/Echo drivers, and can be freely changed during your measurement.

Fourier Step Sum(FSS) technique, which sums ZF-FFT results by changing the center frequency, is employed. The internal memory for the summation is in time domain, which enables you to select window function any time.

“Activate Freq. Sweep” first clears the stored result and start changing a frequency of the SG from “Center Freq.” - “Span”/2 by every “Step”.

“Auto Phase” choses the phase that can maximize the overall real part. However, usually the absolute part should be relevant.


You have to consider “Band Width” for FSS before starting measurement. But you can choose it by multiplying “50%”, “100%”, “200%” anytime.

“LC tuning” determines a method for tuning during the sweep. “As is” does nothing. “Auto Tune” employs an auto-tuner driver and tunes the circuit every “Step”. “Await” turns off the pulser every “Step”, asks you, and waits until you turn on the pulse again after your manual tuning.

## NMR Spectrum Measurement (Field sweep)

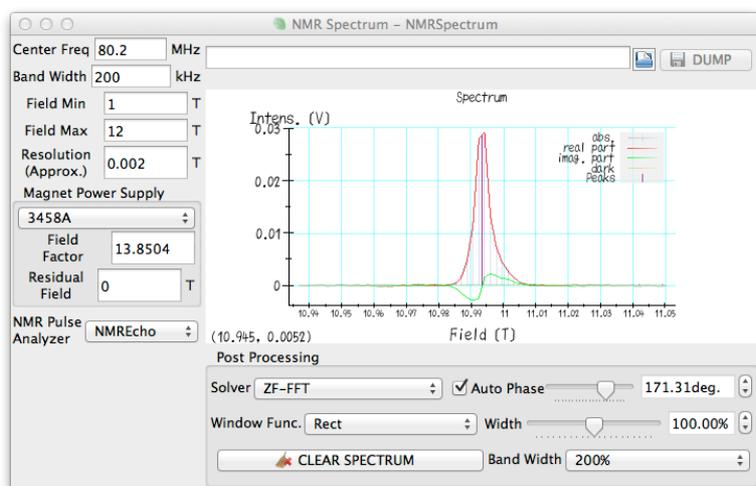This driver is an analyser which measures frequency-swept spectrum using waves in FID/Echo driver. Values of the magnetic field is taken from a magnet power supply driver or DMM driver. Configurations for window function are not shared with FID/Echo drivers, and can be freely changed during your measurement.

Fourier Step Sum(FSS) technique, which sums ZF-FFT results by changing the center frequency, is employed. The internal memory for the summation is in time domain, which enables you to select window function any time.

See also the frequency sweep driver.

Internally, the center of the frequency is calculated as  $-\log(H) \times \text{“Center Freq”}$  during FSS. It is justified as long as the resonance frequency is proportional to the field, or the case “Center Freq”  $\gg 0.001 \times \text{“Band Width”}$ .


The actual field is calculated as (raw value)  $\times \text{“Field Factor”} + \text{“Residual Field”}$ .

You must sweep the magnetic field externally somehow.

“Resolution (Approx.)” specifies a typical resolution of the horizontal axis. Change of this entity may clear the result.

# NMR LC autotuner

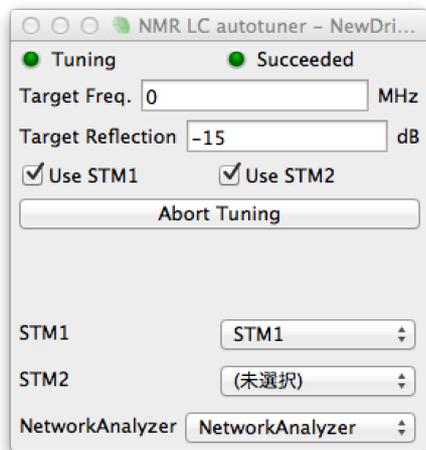This driver is an analyzer, which tunes the LC circuit by monitoring a vector network analyzer driver. Before and after the tuning, it inverts the bits of "AUX" in the motor drivers, which can be used to control an RF relay.

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